

Operating instructions for LABC-Cold-Air-Dryer for TLC records



The LABC-Cold-Air-Dryer serves for the reproduceable dry of TLC records and foils.

Operation:

1. Device drawer forwards pulled out.
2. Record insert concentric and close the drawer again.
3. Net plugs pocket.
4. Device about the green seesaw counter switch on which shines then green.
5. About the black step counter you choose the strength of the airflow.

Step 1: two ventilators

Step 2: four ventilators

Aerial throughput per ventilator (free-blowing): 59 m³/H

Exchanging the filter mat (see also device):

1. Net plugs move.
2. Knurled nut in the back unscrew.
3. Upper case part approx. 20 cm to the back move and take off upwards.
4. The filter mat exchange.
5. Case upper top again from the back put on, up to push all the way to the stop and screw together.
6. Net plugs pocket.

The change interval is directed after the duration of the application time and the lab circumstances. Yearly change is recommended.

All electric components are CE-compliant.

Technical dates: No.: 711862

Operating voltage: 230 V / 50 hertz, 11 watts

For record size: max. 20 x 20 cm

Measures (B x H x T): approx. 23 x 9 x 27.5 cm

Cable length: approx. 2 m